



THS/PH21[®] SERIES

Quality Control at its finest



PHARMACEUTICAL METAL DETECTION SYSTEMS

FEATURES

- Superior and Consistent Detection
- Enhanced Product Throughput Rate
- Preventive Maintenance Not Required
- Maximum Configurability / Integration
- FDA 21 CFR Part 11 Compliant
 - Data Security
 - Data Integrity
 - Data Traceability
- Design and Construction compliant with GMP criteria
- AISI 316L stainless steel construction



- | | | | |
|--|--|--|--|
| <input checked="" type="checkbox"/> TABLETS | <input checked="" type="checkbox"/> GEL CAPS | <input checked="" type="checkbox"/> GRANULAR | |
| <input checked="" type="checkbox"/> CAPSULES | <input checked="" type="checkbox"/> POWDERS | <input checked="" type="checkbox"/> LIQUIDS | <input checked="" type="checkbox"/> SYRUPS |



www.ceia.net



phone +39 0575 4181 • fax +39 0575 418296 • qa-detectors@ceia-spa.com

CEIA THS/PH21 Metal Detection Systems offer detection, construction quality and reliability characteristics that make them the most suitable and effective solution to automatic elimination of metal contaminants.



Validation Solutions Available: Calibration Standards • Validation Manuals
 • Comprehensive Service & Support • Compliance Tools • On-line Production Data Validation

FDA 21 PART 11 COMPLIANT

The THS/PH21 Metal Detector Series is a high-sensitivity, high-precision measuring instrument. The response from the device is used to directly command ejection of the contaminated product and to verify that it has in fact been eliminated.

FDA Code Title 21 Part 11 prescribes rigorous criteria for access to programming and computer data protection, which have been fully adopted in the CEIA THS/PH21 Series firmware. User access is only possible via a User Name and Password login under the control of the System Administrator, who programs and manages individual user profiles and user functions. As a result, the requirements for security, integrity and traceability are satisfied.

All events performed by the THS/PH21 Metal Detector are recorded in a data log to comply with FDA Part 11 protocol and assist in root cause analysis. The events include detection, ejection, program operations, passwords and periodic functional test phases using standard test samples.

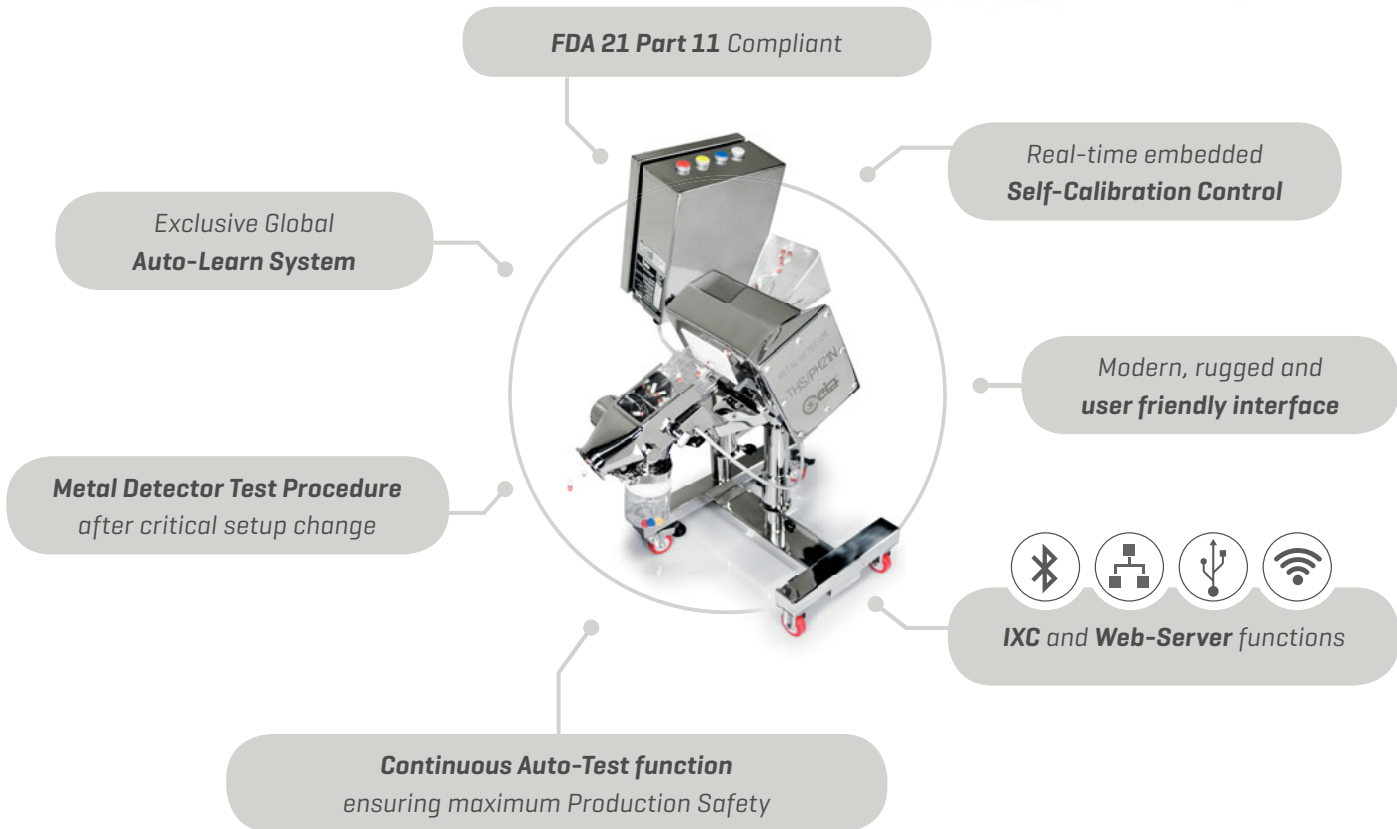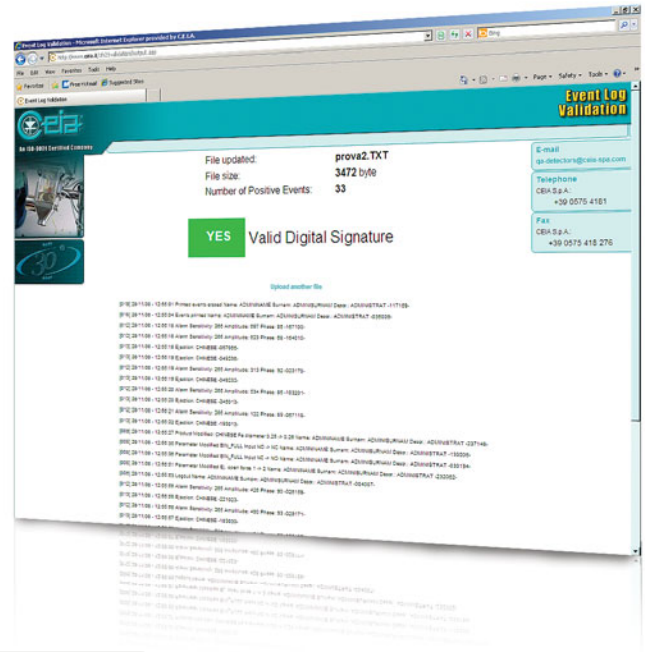

CEIA Metal Detectors are specially engineered so that **no manual calibration or tuning is needed.**

CEIA Metal Detectors have an **internal Auto-diagnosis** for automatic balancing of the antenna with continuous embedded **Self-Calibration control.**

REAL-TIME EMBEDDED SELF-CALIBRATION CONTROL

The Self-Calibration control allows **maximum repeatability and performance consistency over time and with environmental changes.**

By means of specific signals sent to the transmission and reception chain of the Metal Detector, constant monitoring of the detection characteristics is carried out, with consequent constant compensation for any variations caused by environmental factors. The result is system **stability and constancy in detection performance and product effect neutralization.**

CONTINUOUS AUTO-TEST FUNCTION ENSURING MAXIMUM PRODUCTION SAFETY

Special electronic stimuli are sent to the transmission and reception chain of the THS/PH21 Series Metal Detector, causing variations in the detection signals which provide checks on detection characteristics.

These variations are compared with the reference values stored in the memory of the metal detector at the time of the factory calibration test. This produces an automatic certified check of the system's sensitivity.




CEIA offers samples for quality assurance testing certified in composition, size, and electromagnetic response.

EXCLUSIVE GLOBAL AUTO-LEARN SYSTEM

The Auto-Learn System provides **simultaneous maximum sensitivity to all metals starting from a single learning transit.** The THS/PH21 Series employs an exclusive Auto-Learn system for pharmaceutical products, which allows optimization of the detection sensitivity to all metals with the maximum speed and precision, equivalent to hundreds of conventional learning transits. This results in levels of precision and efficiency never before obtained.

NEW DIAGNOSTIC TOOLS

The THS/PH21 Series introduces **new tools for the installation and maintenance technician** to measure the environmental compatibility of the metal detector. These are functions that would otherwise have required the intervention of external instrumentation, now implemented in the microcomputer control of the device itself. The measurements include general mechanical and electromagnetic environmental compatibility.

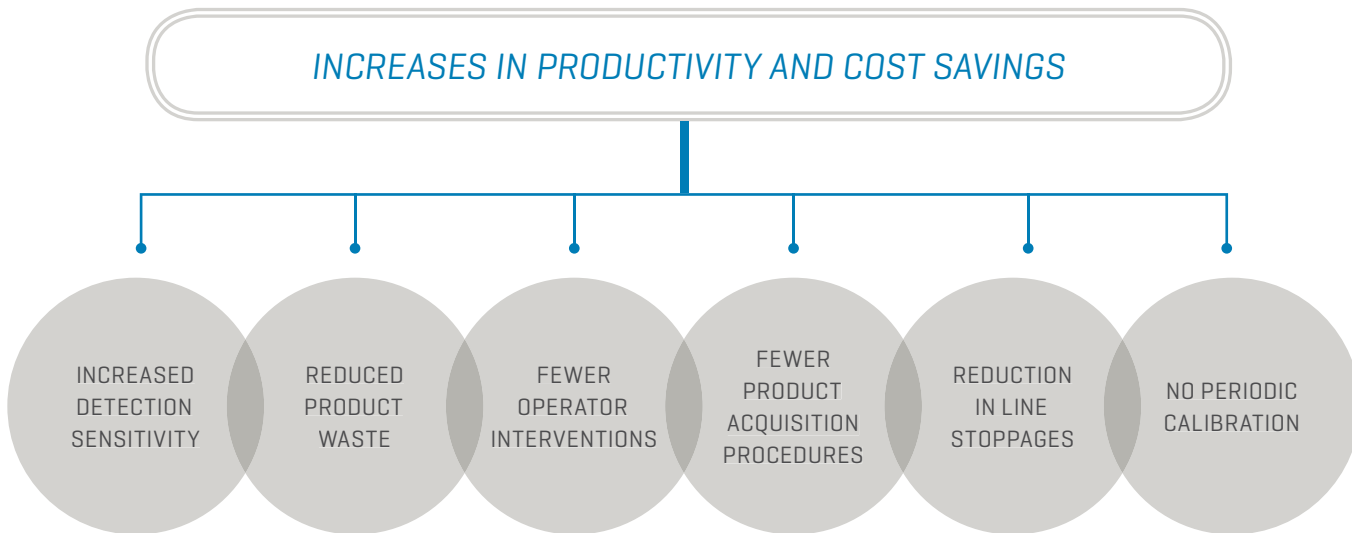


HIGH PERFORMANCE PHARMACEUTICAL METAL DETECTION SYSTEMS



THS/PH21N


ULTRA HIGH SENSITIVITY METAL DETECTOR

- Ultra high sensitivity to all magnetic and non-magnetic metals, including stainless steel
- AISI 316L stainless steel construction
- Control Power Box available according to UL 508A and CSA-C22.2 No. 14-05 
- 500 product profiles, selectable by local programming or network software
- 10,000 storable events
- High-contrast graphic display
- Local programming: 16 stainless steel keys, 3 with double function
- Quick Access key for fast programming of user parameters
- Antistatic, Jamming Sensor and Bin full Kits [available on demand]



THS/PH21E

HIGH SENSITIVITY METAL DETECTOR

- High sensitivity to all magnetic and non-magnetic metals, including stainless steel
- AISI 316L stainless steel construction
- Control Power Box available according to UL 508A and CSA-C22.2 No. 14-05 
- 250 product profiles, selectable by local programming or network software
- 1,000 storable events
- 4 x 20 character alphanumeric display
- Local programming: 4 stainless steel keys, 3 with double function
- Antistatic, Jamming Sensor and Bin full Kits [available on demand]

Due to many cross-platform studies made by CEIA engineers, the THS Metal Detector Series 21 is HIGHLY IMMUNE to common radio frequencies and difficult environments typically found in industrial installations.



THS/PH21N-DT • THS/PH21E-DT

DUST TIGHT METAL DETECTION SYSTEM

Dust tight transit pipe and ejection system prevents dispersion of the product

- Transit tunnel and ejection system completely sealed
- Easy integration with tablet press and deduster from the leading manufacturers
- Rapid maintenance does not require the employment of tools to assemble and disassemble the components
- Antistatic, Jamming Sensor and Bin full Kits [available on demand]



THS/PH21N-WIP • THS/PH21E-WIP

WASH-IN-PLACE METAL DETECTION SYSTEM

A special built-in washing system allows complete cleaning of the conduits after completion of each production batch

- Transit tunnel and ejection system completely sealed
- Fitted with nozzles for washing water injection and decontamination
- Easy integration with tablet press and deduster from the leading manufacturers
- Antistatic, Jamming Sensor and Bin full Kits [available on demand]




SPECIFIC INSPECTION AGAINST TINY METAL PARTICLE CONTAMINATION



THS/PH21N-D25


ULTRA HIGH SENSITIVITY METAL DETECTOR

- Unsurpassed Detection
- Best Product Quality
- Round aperture
- Applications:
 - ✓ R&D analysis
 - ✓ Rework confirmation
 - ✓ Root cause analysis assistance
 - ✓ New product specification

PRODUCT		VERIFIED THROUGHPUT [items / h]
		THS/PH21N-D25
	Diameter: Ø 8,5 mm Thickness: 4,5 mm	30,000

EXAMPLES OF THROUGHPUT

- Tilt Angle: 35°
- Max Throughput Rates are related to the shape and to the size of the tablets

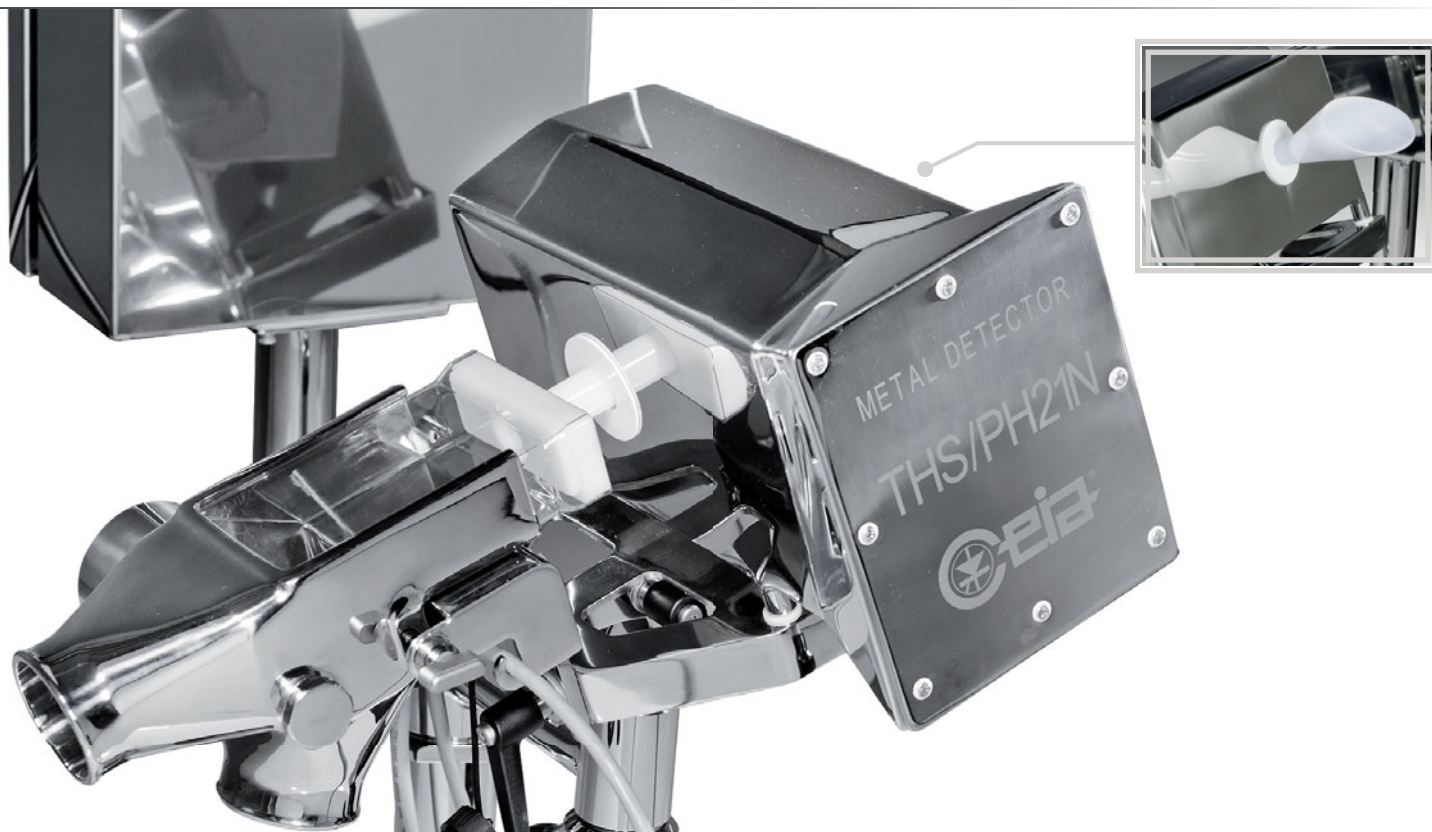

INTEGRATED SYSTEMS FOR GRANULAR, LIQUID AND POWDER PRODUCTS



THS/PH21N-FFV • THS/PH21E-FFV

INTEGRATED SYSTEM FOR **GRANULAR** AND **POWDER PRODUCTS**

The ejection system is fitted with an innovative deflector that features quick response, precise ejection of the contaminated product and high production flow rate.

- Ultra High Sensitivity to all metals
- Very compact size
- AISI 316L stainless steel construction
- Control Power Box available according to UL 508A and CSA-C22.2 No. 14-05
- ATEX ZONE 21 version [available on demand]



THS/FFV 21 SERIES

LARGE APERTURE INTEGRATED SYSTEM FOR **GRANULAR** AND **POWDER PRODUCTS**

- Wide range of standard pipe sizes available to suit all applications
- Enhanced product throughput rate
- AISI 316L stainless steel construction
- Compact size
- Fast and efficient: achieve fast ejection speeds and accuracy, with the minimum amount of rejected product
- Pipe in antistatic material, certified for contact with food products
- ATEX version [available on demand]



THS/PLV 21 SERIES



INTEGRATED SYSTEM FOR **LIQUID** AND **VISCOUS PRODUCTS**

- Wide range of standard pipe sizes available to suit all applications
- Full compliance with HACCP criteria
- AISI 316L stainless steel construction
- Fast acting reject device for a limited quantity of rejected product
- Easy to clean and inspect

THS/PH21® MECHANICAL DATA



THS/PH21N • THS/PH21E models

MODELS	APERTURE SIZE [DW x DH]
THS/PH21N THS/PH21E	90 x 25
	100 x 40
THS/PH21N-DT THS/PH21E-DT	90 x 25
	100 x 40
THS/PH21N-WIP THS/PH21E-WIP	90 x 25
	100 x 40

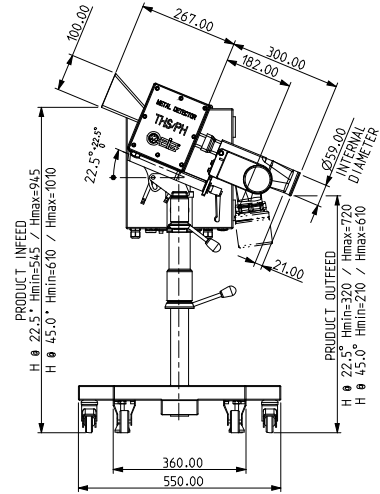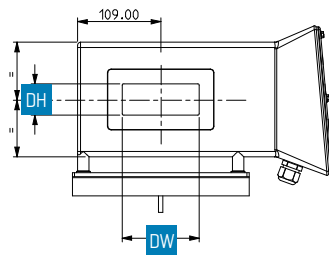


THS/PH21N-D25 model

MODEL	APERTURE SIZE
THS/PH21N-D25	Ø 25

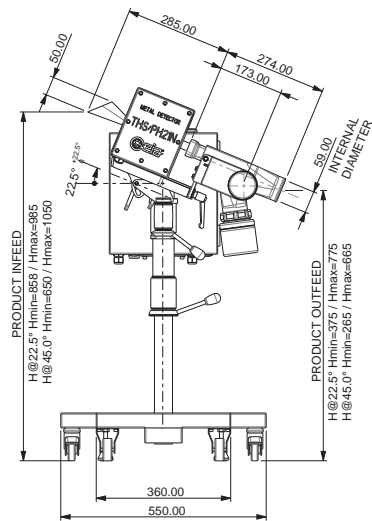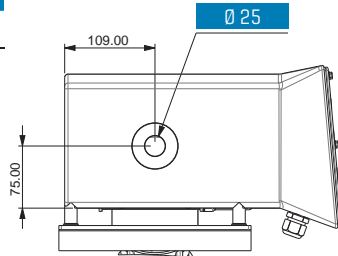

THS/PH21N-FFV • THS/PH21E-FFV models

MODELS	APERTURE SIZE	A	B
THS/PH21N THS/PH21E	100 x 40	90	30
	90 x 25	78	13

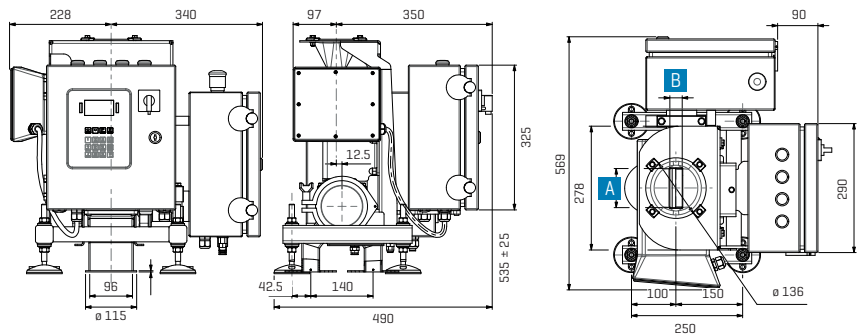


Dimensions in mm

THS 21 METAL DETECTORS SERIES WITH MULTI-SPECTRUM TECHNOLOGY FOR CONVEYOR INSPECTION SYSTEMS



THS/FBB • THS/MBB

FLAT AND MODULAR CONVEYOR INSPECTION SYSTEMS

- High Sensitivity to all metals
- Integrated controls for line speed and rejection
- High immunity to environmental interference
- Automatic learning & tracking of product effect
- Easy to maintain and inspect
- Digitally adjustable belt speed
- Control Power Box available according to UL 508A and CSA-C22.2 No. 14-05 
- Applications:
 - Flacons
 - Bottles
 - Packaging
 - Loose product



BASIC CONFIGURATION

optical signalling device		acoustic signalling device	
belt stop in case of alarm		synchronization photocell	
variable speed		complete reversibility of the accessory assembly and of the Metal Detector	
conveyor belt dimensions (LxWxH)		1800 mm x 85 mm x 875 mm	1800 mm x 150 mm x 875 mm
type of stand	feet	●	●
	castors	○	○
flap reject		○	○
ejection confirmation/bin full kit		○	○
insufficient air pressure control kit		○	○
side panels for product containment		●	●
test due signalling device [blue]		○	○
● STANDARD ○ ACCESSORIES / OPTIONALS			



THS/PH21® ADVANTAGES AND PERFORMANCES



MODERN, RUGGED AND USER FRIENDLY INTERFACE

- INDUSTRIAL RATE DESIGN
- RAPID DATA ENTRY WITH EXTENDED KEYBOARD
- EASY TO READ, HIGH-CONTRAST GRAPHIC DISPLAY
- RUGGED, ANTIVANDALIC STAINLESS STEEL KEYBOARD

Organic graph display with extremely high contrast [3000:1] and a viewing angle up to 180°, and an extended alphanumeric keypad for quick keystrokes.

- PROTECTION DEGREE:
 - Power Supply Box: IP65 [IEC60529]
 - Antenna & Ejection System: IP67
- 40 USERS AVAILABLE [INCLUDING ADMINISTRATOR]
- QUICK ACCESS FUNCTION


A new function, activated simply by pressing a dedicated key, allows the most frequently-used functions to be directly recalled. These functions are programmable by the operator.



EJECTION SYSTEM SPECIFICATIONS

- Monolithic AISI 316L Stainless Steel Body
- Advanced mechanical and electronic solutions guarantee the integrity of the inspected product
- Extremely quick response time provides full safety in the ejection of contaminated product
- Double confirmation system carefully monitors and controls the complete opening and closing of the ejection flap
- Bin full sensor kit [on request]

- 1 Metal Detector Ejection System
- 2 Metal Detector Sampling Ejection and Safe Test
- 3 Product Sampling
- 4 Contaminated Product

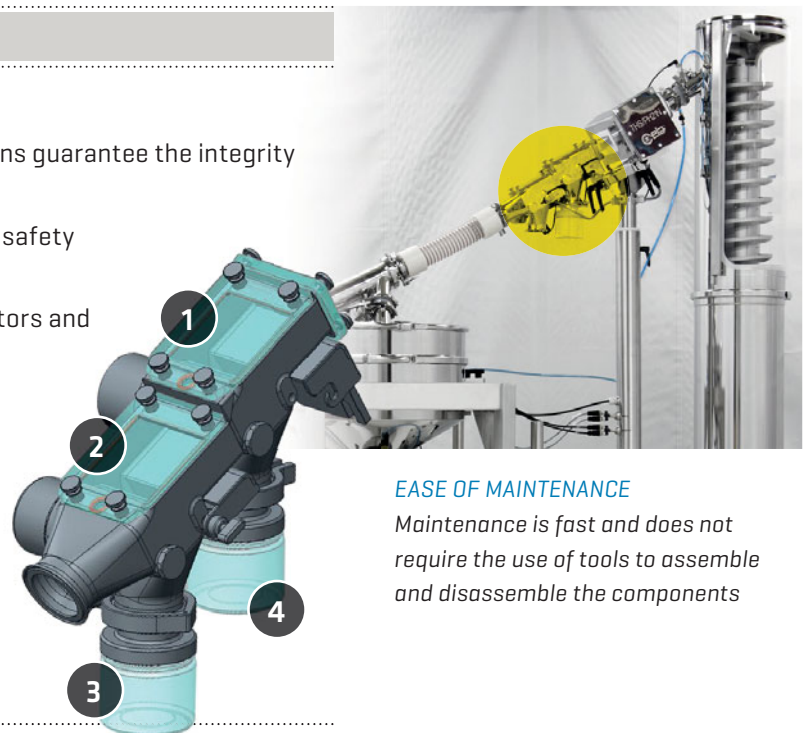

EASE OF MAINTENANCE

Maintenance is fast and does not require the use of tools to assemble and disassemble the components


SOPHISTICATED INTERFACE CAPABILITIES

BLUETOOTH® CONNECTIVITY

In the THS/PH21 Series, local connection to the maintenance technician’s computer no longer requires physical access to the interior of the detector or the use of unwieldy connection cables. The Bluetooth connection can be used for programming, monitoring of the signals via the CEIA MD-Scope program and the transfer of the data contained in the Metal Detector’s events memory.

NETWORK COMMUNICATION CAPABILITIES

The Metal Detector can be linked to an Ethernet network (optional IXC module required). In association with the THS Production Plus Software it enables remote management of production, collection of all technical and operational events, generation of statistical and traceability reports in compliance with FDA 21 CFR part 11 requirements.

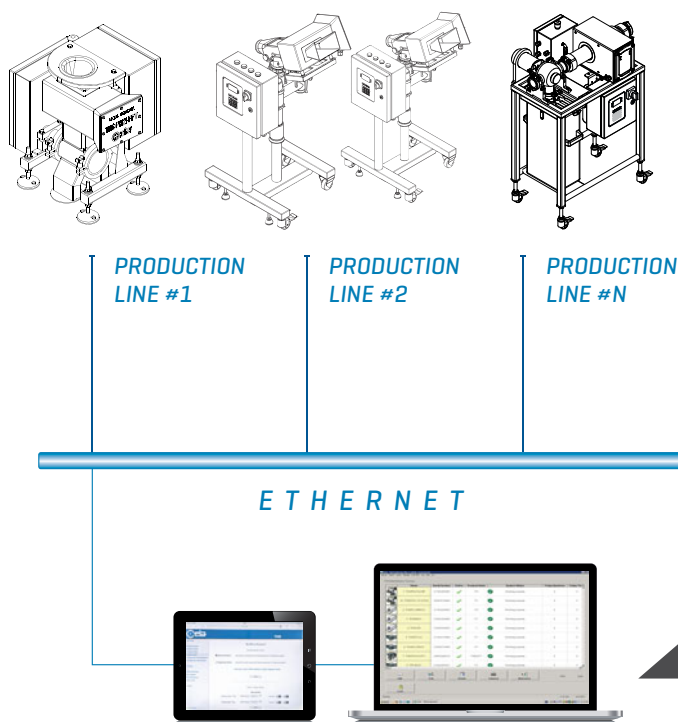

THS PRODUCTION PLUS

MD-SCOPE

Software Diagnostic Package complete with connector cable and hardware key:

- Remote programming
- Signal input/output
- Alarm threshold and amplitude of the signal
- Oscilloscope functionalities

THS PRODUCTION PLUS

- Connects and Acquires Data from Multiple THS Detectors via Ethernet LAN
- Provides Connection to an External Database and Definition of a Programmable Block of SQL Instructions for each Metal Detector Event
- Enables Data Base Management and Back-Up of Metal Detector events
- Report Data Exportable in HTML and CSV format
- Provides User Authentication and Manages Electronic Signatures and Records
- Allows Data Integrity and Operator Auditing
- FDA Requirements Compliant, Title 21 of the Code of Federal Regulations [CFR]
- E-mail sender

Name	Serial Number	Online	Product Name	System Status	Serial Number	Today Tests	Last S
1 THSPH21N-WIP	21100208062	✓	P2	Working properly	0	28/20	
3 THSMS21	2100024065	✓	DEFAULT	Working properly	0	28/20	
4 THS21E	21000238032	✓	Y	Working properly	0	28/20	
5 THSFFV21	21100213069	✓	A	Working properly	0	0	28/20
6 THSPLVMS21	21000218048	✓	P1	Working properly	0	0	28/20
7 THSPH21N-FFV	20900246015	✓	TABLET1	Working properly	0	0	28/20
8 TPLMS21	21000205021	✓	P1	Working properly	0	0	28/20
Total:							

IXC AND WEB-SERVER FUNCTIONS



IXC MODULE

- Gateway to THS Production Net Management System
- Integrated Web-server
- Dual 100base-T Ethernet
- Dual Full Speed USB interface
- Wi-Fi 802.11b/g
- 100,000 events internal data logger
- Additional web-server functionality

LOCAL CONNECTIVITY




ADVANCED CONNECTIVITY



IXC
MODULE

NETWORK

WINDOWS BASED SERVER WITH NETWORK SEARCH ENGINE

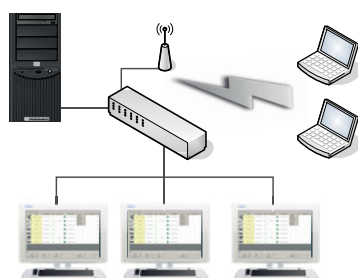

WIRED NETWORK CLIENTS

WIRELESS NETWORK CLIENTS

WEB-SERVER INTERFACE

BUILT-IN MD SCOPE

The MD Scope functions (including oscilloscopes) are available to every user connected to the Metal Detector via Ethernet or WiFi.

STATUS

Replicates every message displayed by the THS (updated each 5 s).

PRODUCTS

Displays the list of products stored into the THS memory, the selected product is shown in green. This page allows the product to be changed.

PRODUCTS EXPORT

Allows selection of all or part of the products stored in the THS memory in order to export them.

PRODUCTS IMPORT

Allows importation of all or part of the products exported.

BUFFER EXPORT

Allows exportation of events occurred in a selected time frame.

CONFIGURATION BACKUP/RESTORE

Allows backup or restoring of the connected system configuration.

CONFIGURATION EXPORT/IMPORT

Allows the configuration to be exported or imported from other Metal Detectors of the same model, installed on the same system.



CEIA COSTRUZIONI ELETTRONICHE AUTOMATISMI INDUSTRIALI

Zona Industriale 54/G, 52041 Vicinaggio - Arezzo [ITALY]

Tel.: +39 0575 4181 Fax: +39 0575 418296 E-mail: pharma@ceia-spa.com

CÔNG TY TNHH CÔNG NGHỆ HOÀNG LONG

468 Lê Hồng Phong, P.1, Q.10, Tp. HCM

Hotline: 0907.686.889

info@congnghehoanglong.com

www.congnghehoanglong.com